
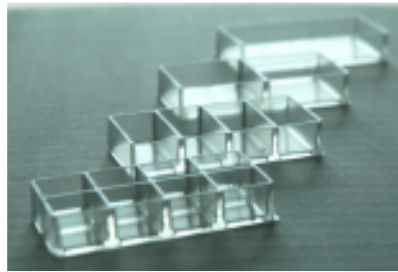
Imaging Chamber / Imaging Chamber CG

Instructions for handling and chamber
disassembly

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
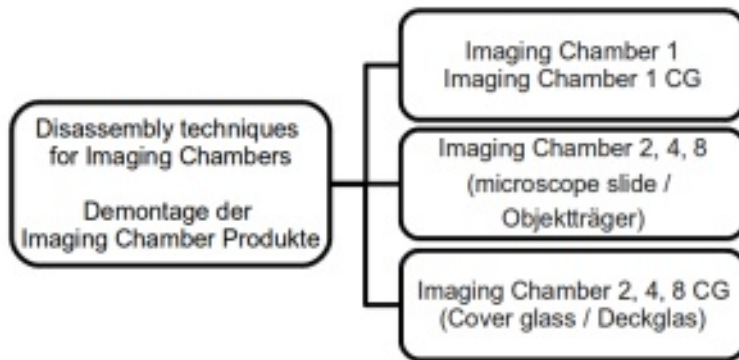
All Imaging Chamber and Imaging Chamber CG products are designed to allow disassembly of the chamber body and the glass bottom at the end of experiments / investigations. Due to their physical characteristics different strategies for microscope slide bottoms and the cover glass bottoms have to be considered. Also the chamber design (one well versus multiwell) determines the optimal technique for disassembly.

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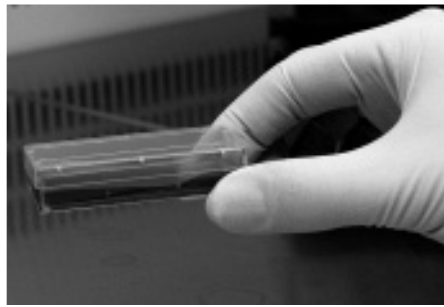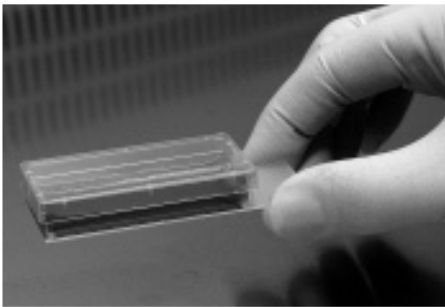
Handling / Handhabung

During transport we recommend to hold the chambers where they are supported by a wall or to grip the microscopic slide (Imaging Chamber products).

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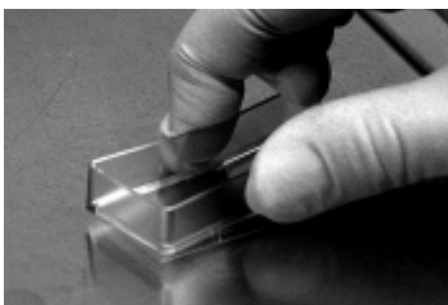Disassembly / Demontage Imaging Chamber 1 und Imaging Chamber 1 CG


To disassemble the glass bottom from Imaging Chamber 1 or Imaging Chamber 1 CG you only need to compress the two sidewalls of the chamber. The shear force detaches the glass from the chamber.

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



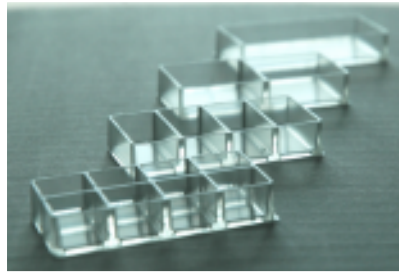
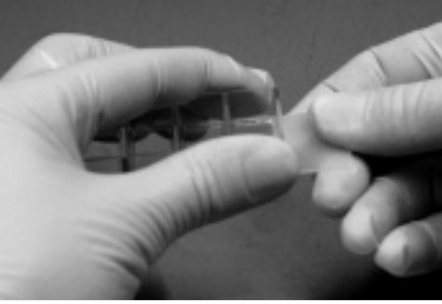
To disassemble the microscope slide from Imaging Chamber 2, 4 or 8 you have to hold the slide at the outer frosted end close to the chamber while lifting the chamber beginning from this end with the other hand.

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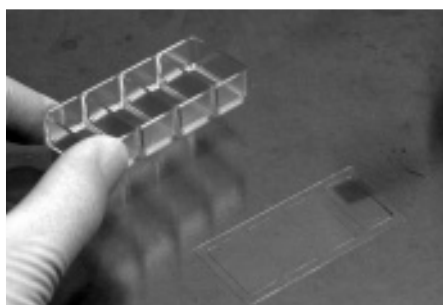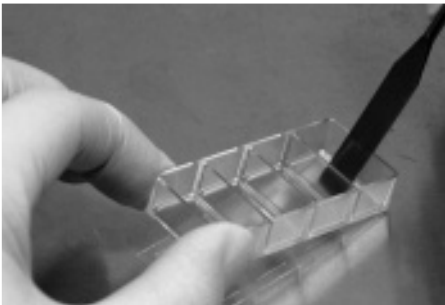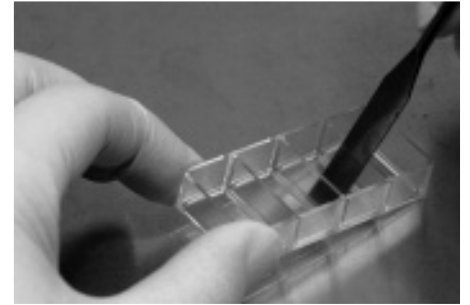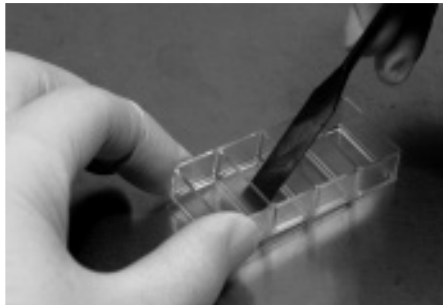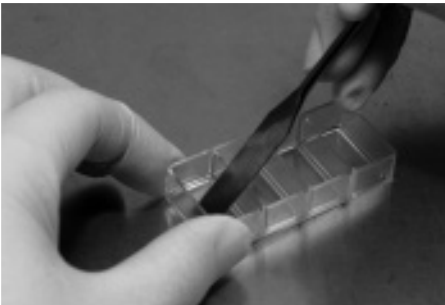Disassembly / Demontage Imaging Chamber 2, 4 und 8 CG Cover Glass / Deckglas

To disassemble the cover glass from Imaging Chamber 2, 4 or 8 CG you will need a tool like a spatula. Place the imaging Chamber on a clean and flat ground. Lift the Chamber approximately 1 mm at one end of the Chamber from the ground. Gently press the cover glass with the spatula away from the chamber body close to the inner Chamber walls. Proceed well by well until the glass comes off.

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Video Tutorials can be found at www.zell-kontakt.de

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